# **In**Spec<sup>™</sup> Data Sheet

### **Cleanroom Contamination Control**



## Manual Detergent Pre-Diluted Solution

An alkaline detergent and hard surface cleaner for use in the biopharmaceutical and cleanroom industries.



### GENERAL GUIDANCE

#### USE INSTRUCTONS: Do not dilute.

InSpec N10 is designed for cleaning hard surfaces. Remove any outer packaging and spray onto the surface or pour into a suitable bucket. Apply the solution onto the surface to be cleaned. Ensure good cover of the surface with the solution. Clean the surface with a mop or wipe. Rinse or squeegee away residue as required.

#### TECHNICAL DATA:

N10 is an alkaline solution (pH 12) containing 4% actives. The product is a blend of alkaline detergent, surfactant and sequestering agents designed to assist the removal of process residue and general solling.

#### STORAGE:

Store in original packaging, away from sunlight and extremes of temperature. Full guidance on the handling and disposal of this product is provided in the Material Safety Data Sheet (MSDS).

#### SAFE HANDLING:

Always wear gloves and goggles or face protection. Always read the label and the MSDS before use.

#### MATERIAL COMPATIBILITY:

Application of solutions, when used as directed, will not affect materials normally encountered in the cleanroom. See compatibility information in the Technical File.





#### BENEFITS

- Designed for cleaning floors, walls, ceiling and other hard surfaces.
- The Dycem recommended cleaner of Dycem flooring.
- Rapid action alkaline detergent.
- Suitable for manual cleaning.
- Available in convenient ready to use formats.
- Designed for the removal of general process derived residue and soiling.
- A useful solution for removing residue accumulated by the regular use of rotational biocides.
- Defined 24 month shelf life.
- InSpec N10 trigger spray has a validated 3 month in-use shelf life.

#### **TECHNICAL USER INFORMATION**

InSpec N10 is a pre-diluted detergent.

Prediluted formats use WFI quality water.

InSpec N10 is manufactured to GMP in an ISO 6 cleanroom. The product is filled and capped in an ISO class 5 environment.

The solution is filtered through a 0.2 micron filter into the final packaging.

All InSpec products are suppled with key certification such as sterility and conformity.

#### InSpec N10 PRODUCTS

PRODUCT	PACK SIZE	CARTON SIZE	PRODUCT CODE	STERILITY
InSpec™N10 1L Trigger	1L trigger spray	6 X 1L	N10WFI1CC-1LS	Sterile
InSpec™N10 1L Trigger	1L trigger spray	6 X 1L	ICN10WFI-1NS	Non-Sterile
InSpec™N10 5L Screw Cap	5L screw cap	2 X 5L	N10WFI1-5LS	Sterile
InSpec™N10 5L Screw Cap	5L screw cap	2 X 5L	ICN10WFI1-5NS	Non-Sterile
InSpec™N10 Concentrate	100ml concentrate	50 X 100ml	N10WFI1-100MLS	5 Sterile
InSpec <sup>™</sup> N10 Burstable Wipes	15 wipes/pouch	20 pouches	N10WFI60BP15S	Sterile

Discovery 2 • NETPark • 2 William Armstrong Way • Sedgefield • County Durham • TS21 3FD • UK UK Tel: 0845 073 9529 • International Tel: +44 (0) 1527 830 942 Web: www.redditchmedical.com • Email: sales@redditchmedical.com